


C3000 Replacement Syringes

Cat#	10070	10071	10072	10073	10074	10075	10076
Volume	50µL	100µL	250µL	500µL	1mL	2.5mL	5mL
Length	3cm	3cm	3cm	3cm	3cm	3cm	3cm
Fitting Interface	¼ - 28	¼ - 28	¼ - 28	¼ - 28	¼ - 28	¼ - 28	¼ - 28
Materials	UHMWPE seal, Borosilicate Glass	UHMWPE seal, Borosilicate Glass	UHMWPE seal, Borosilicate Glass	UHMWPE seal, Borosilicate Glass	UHMWPE seal, Borosilicate Glass	UHMWPE seal, Borosilicate Glass	UHMWPE seal, Borosilicate Glass

Cat#	10080	10081	10082	10083	10084	10085	10086
Volume	50µL	100µL	250µL	500µL	1mL	2.5mL	5mL
Length	3cm	3cm	3cm	3cm	3cm	3cm	3cm
Fitting Interface	¼ - 28	¼ - 28	¼ - 28	¼ - 28	¼ - 28	¼ - 28	¼ - 28
Materials	PTFE seal, Borosilicate Glass	PTFE seal, Borosilicate Glass	PTFE seal, Borosilicate Glass	PTFE seal, Borosilicate Glass	PTFE seal, Borosilicate Glass	PTFE seal, Borosilicate Glass	PTFE seal, Borosilicate Glass

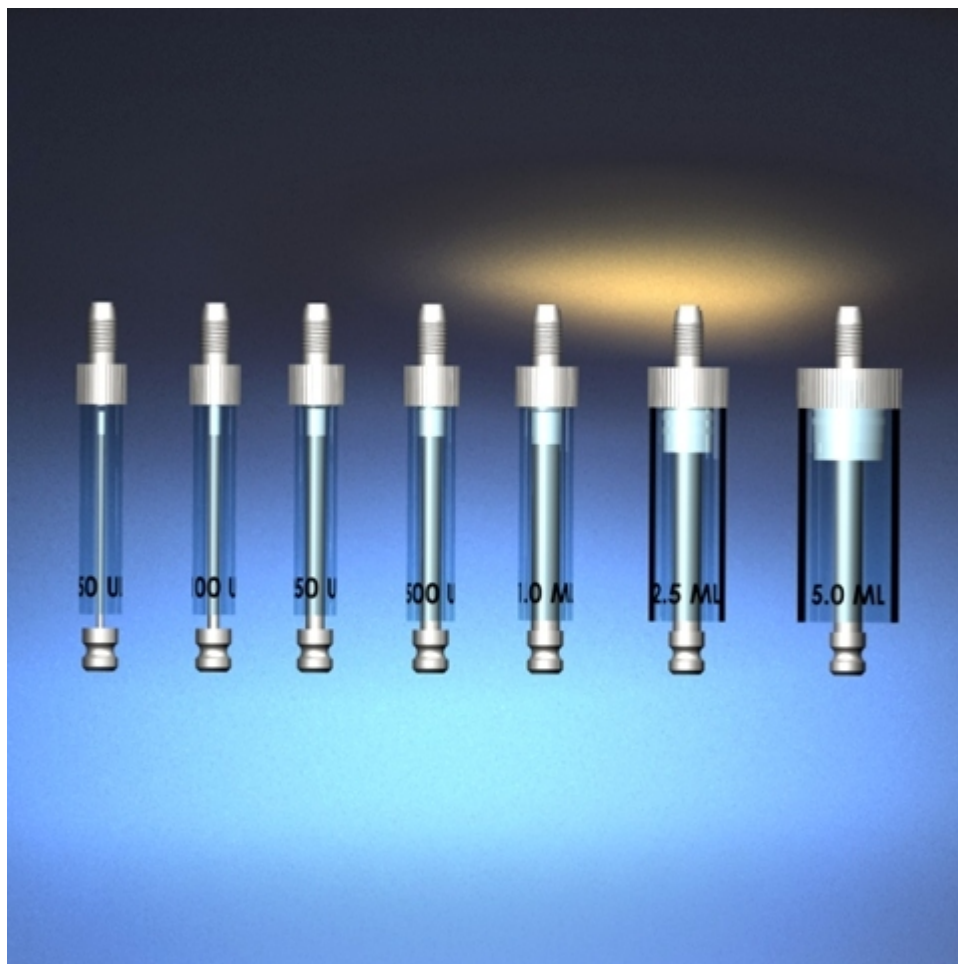

SYRINGE REPLACEMENT

To replace the syringe, prime all fluid from the syringe, and follow the steps below:

1. To replace the syringe, lower the plunger driver. If the power is on, this can be done by sending the A3000R command. If the pump is powered off, manually lower the plunger by pushing firmly down on the syringe plunger holder.
2. Turn the plunger lock screw counter clockwise 3 turns.
3. Turn the syringe counter clockwise and remove from the valve.

To install the syringe follow the steps below:

4. Screw the threaded portion of the syringe clockwise into the valve
5. Fully seat the plunger into the syringe plunger holder.
6. Securely tighten (clockwise) the plunger lock screw.
7. Initialize the pump.

